

DeliverX™ siRNA Transfection Evaluation Kit 0.12 mL

Product Insert

Cat. No. DX0001

**About
DeliverX siRNA
Evaluation Kit**

The DeliverX siRNA Transfection Evaluation Kit includes all of the components to enable a new user to evaluate our DeliverX siRNA solution for efficiency of siRNA transfection and simple assay workflow for a cell type of interest. The evaluation kit includes the DeliverX siRNA Transfection Kit and two proficiency controls, FAM-labeled siRNA Control and Human GAPDH siRNA Control, that enable the user to optimize transfections under controlled conditions.

**About
DeliverX siRNA
Transfection Kit**

Panomics' DeliverX siRNA Transfection Kit is suitable for the transfection of most cell types. The kit contains the reagents required to efficiently transfect siRNAs into most cell types with minimal damage and good reproducibility when optimization guidelines provided in the User Manual are followed.

**About
FAM-Labeled
siRNA Control**

Panomics' FAM-labeled siRNA Control enables visualization (1–4 hours post transfection) of transfection efficiency. This validated siRNA delivery control is a 21-mer, double-stranded, deoxyoligonucleotide.

**About
Human GAPDH
siRNA Control**

Panomics' Human GAPDH siRNA Control is a validated potent siRNA for functional evaluation of transfection efficiency in human cell types.

**Safety Warnings
and Precautions**

All chemicals should be considered potentially hazardous. We recommend that this product and its components be handled by those trained in laboratory techniques and be used according to the principles of good laboratory practice.

For research use only. Not for use in diagnosis of disease in human or animals.

**Contents and
Storage**

Component	Volume	Storage
DeliverX siRNA Transfection Reagent	0.12 mL	-20 °C
FAM-Labeled siRNA Control	0.12 mL	-20 °C/Dark
Human GAPDH siRNA Control	0.06 mL	-20 °C
DeliverX siRNA Buffer-1	4 mL	15–30 °C
DeliverX siRNA Buffer-2	4 mL	15–30 °C

If stored properly, product shelf life is 6 months from the date received.

**Contacting
Panomics**

For technical questions, contact our technical support group by telephone at 1-877-726-6642 option 3 or by email at techsupport@panomics.com (US and Canada) techsupport_europe@panomics.com (Europe), or visit our website www.panomics.com for an updated list of FAQs and product support literature.

DeliverX is a trademark of Panomics, Inc.

DeliverX and DeliverX Plus transfection reagents are manufactured and distributed by Panomics, under license from CNRS (France).

© 2006 Panomics, Inc. All rights reserved.

Panomics Inc., 6519 Dumbarton Circle, Fremont, CA 94555, Tel. (877) 726-6642
